

Alexa Fluor^{®*} 647 FLUORONANOGOLD^{™†} Conjugates



95 Horseblock Road, Unit 1, Yaphank NY 11980-9710
Tel: (877) 447-6266 (Toll-Free in US) or (631) 205-9490 Fax: (631) 205-9493
Tech Support: (631) 205-9492 tech@nanoprobes.com
www.nanoprobes.com

PRODUCT INFORMATION

ALEXA FLUOR[®] 647 FLUORONANOGOLD[™]-ANTIBODY CONJUGATES

Catalog Numbers & Product Names:	7501	Alexa Fluor [®] 647 FluoroNanogold [™] goat anti-mouse IgG (H+L)
	7502	Alexa Fluor [®] 647 FluoroNanogold [™] Fab' fragment of goat anti-mouse IgG (H+L)
	7503	Alexa Fluor [®] 647 FluoroNanogold [™] goat anti-rabbit IgG (H+L)
	7504	Alexa Fluor [®] 647 FluoroNanogold [™] Fab' fragment of goat anti-rabbit IgG (H+L)
	7505	Alexa Fluor [®] 647 FluoroNanogold [™] rabbit anti-goat IgG (H+L)
	7506	Alexa Fluor [®] 647 FluoroNanogold [™] Fab' fragment of rabbit anti-goat IgG (H+L)
	7507	Alexa Fluor [®] 647 FluoroNanogold [™] goat anti-rat IgG (H+L)
	7508	Alexa Fluor [®] 647 FluoroNanogold [™] Fab' fragment of goat anti-rat IgG (H+L)
	7552	Alexa Fluor [®] 647 FluoroNanogold [™] goat anti-human IgG (H+L)
	7554	Alexa Fluor [®] 647 FluoroNanogold [™] goat anti-guinea pig IgG (H+L)
Appearance:		Pale turquoise liquid
Storage:		Upon receipt store product at 2-8°C. Product is shipped at ambient temperature.
Revision:		1.0 (August 2016)

Nanoprobes offers a new line of Alexa Fluor[®] 647 FluoroNanogold[™] antibody and streptavidin conjugates labeled with both Alexa Fluor[®] 647 dye and the 1.4 nm Nanogold[®] particle. The dual labels are covalently attached to the antibody or streptavidin, enabling both fluorescence and electron microscope observation of the same structure in a single labeling procedure. The Alexa Fluor[®] 647 dye is a bright, far-red-fluorescent dye with excitation ideally suited for the 594 nm or 633 nm laser lines, and is useful in stochastic optical reconstruction microscopy (STORM). It can also be used in two-photon excitation (TPE) microscopy. The presence of the 1.4 nm Nanogold[®] particle allows electron or light microscope detection with and without silver or gold enhancement. The presence of both the fluorescent and the gold labels make correlative microscopy study possible.

* Alexa Fluor is a registered trademark of Molecular Probes, Inc. / Thermo Fisher Scientific, Inc.

† Patented technology.

CONTENTS

Product Information

General Considerations for Immunostaining with Alexa Fluor[®] 647 FLUORONANOGOLD[™] Conjugates

Using Stains with Alexa Fluor[®] 647 FLUORONANOGOLD[™]

Temperature Caution

Thiol Caution

Fluorescence Microscopy with Alexa Fluor[®] 647 FLUORONANOGOLD[™]

Superresolution Microscopy with Alexa Fluor[®] 647 FLUORONANOGOLD[™]

Electron Microscopy Immunolabeling with Alexa Fluor[®] 647 FLUORONANOGOLD[™]


1. Cells in Suspension
2. Thin Sections

Special Considerations for Viewing Alexa Fluor® 647 FLUORONANOGOLD™ in the Electron Microscope
 Silver Enhancement of Alexa Fluor® 647 FLUORONANOGOLD™ for EM
 Gold Enhancement of Alexa Fluor® 647 FLUORONANOGOLD™ for EM
 References

Warning: For research use only. Not recommended or intended for diagnosis of disease in humans or animals. Do not use internally or externally in humans or animals. Non radioactive and non carcinogenic.

PRODUCT INFORMATION

Fab' Conjugate

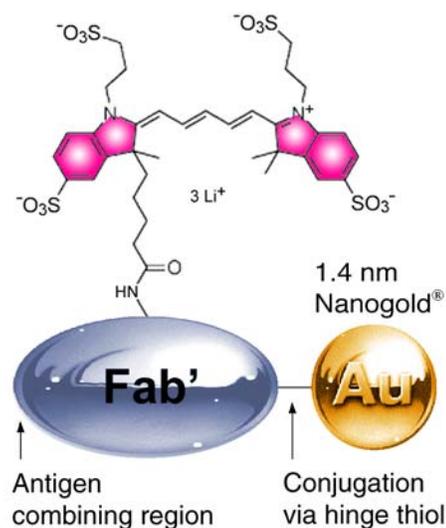

IgG Conjugate

Available only for
 Alexa Fluor® 647
 FluoroNanogold™

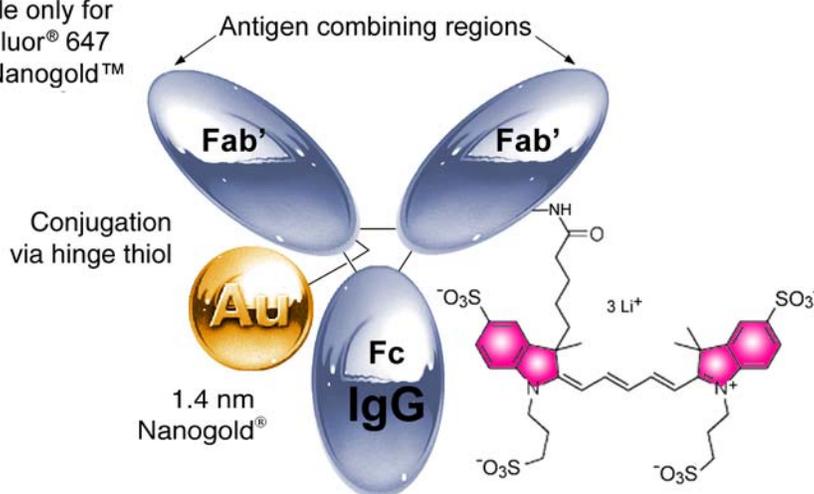


Figure 1: IgG and Fab' covalently conjugated to Alexa Fluor® 647 and NANOGOLD® to give Alexa Fluor® 647 FLUORONANOGOLD™.

Alexa Fluor® 647 FLUORONANOGOLD™ is a unique, dual-purpose probe¹. Alexa Fluor® 647 FLUORONANOGOLD™ conjugates consist of affinity-purified IgG molecules, Fab' fragments or streptavidin conjugated to both Alexa Fluor® 647 dye and the 1.4 nm NANOGOLD® particle.² The structures of Alexa Fluor® 647 FLUORONANOGOLD™ IgG and Fab' conjugates are shown in Figure 1. In the fluorescence microscope, these probes may be used just like conventional fluorescently-labeled antibodies,³ while in the electron microscope they are visualized in exactly the same manner as for NANOGOLD® conjugates⁴. The combination of fluorescence and electron microscopy, which is referred to as correlative microscopy¹³, allows imaging of the same structure in both microscopes. The covalent label linkage is stable indefinitely, and the attachment of NANOGOLD® at a hinge thiol site ensures maximum preservation of native immunoreactivity. These reagents are supplied at a concentration of 0.08 mg/mL of antibody dissolved in 20 mM phosphate buffered saline (150 mM NaCl) at pH 7.4 (PBS), with 0.1 % BSA and 0.05 % sodium azide as preservatives. Alexa Fluor® 647 FLUORONANOGOLD™ conjugates should be stored at 2-8°C. DO NOT FREEZE. PROTECT FROM LIGHT.

GENERAL CONSIDERATIONS FOR IMMUNOSTAINING WITH FLUORONANOGOLD™ REAGENTS

Basically, normal methodologies for each component of the label may be used successfully with Alexa Fluor® 647 FLUORONANOGOLD™ labeling agents. Due to some quenching of fluorescence by the gold particle, slightly higher concentrations of antibody are recommended for incubations. A blocking agent of 5% non-fat dried milk has been found to reduce nonspecific background staining in some cases: this should be used before incubation with probe (in standard wash/blocking steps), and additionally, the Alexa Fluor® 647 FLUORONANOGOLD™ probe can be diluted in a solution also containing 1% non-fat dried milk

before it is applied. The nonspecific staining can also be reduced or eliminated by using a biotinylated antibody as a secondary and a FLUORONANOGOLD™-Streptavidin as a tertiary probe.

PROPERTIES

Alexa Fluor® 647 FLUORONANOGOLD™ contains an extremely uniform 1.4 nm diameter gold particle ($\pm 10\%$), and 3-6 fluorophores per one IgG, or 2-3 fluorophores per one Fab'. The exact degrees of labeling are reported on the specification sheet for each product lot. The absorption maximum of Alexa Fluor® 647 occurs near 650 nm, and the emission maximum is near 668 nm. These values are very similar to those of the Cy5 dye. The Alexa Fluor® 647 dye is readily excited by the 594 nm or 633 nm laser lines.

Alexa Fluor® 647 FLUORONANOGOLD™ contains 1.4 nm NANOGOLD® attached at the antibody hinge region. The small diameter gold particle provides the least quenching of the fluorescent signal from nearby fluorophores, and makes the conjugate the smallest gold immunoprobe commercially available. FLUORONANOGOLD™ will penetrate and reach antigens inaccessible to larger colloidal gold probes.

IgG - Alexa Fluor® 647 FLUORONANOGOLD™ provides a bright fluorescent signal and very low nonspecific staining.

Fab' - Fluor® 647 FLUORONANOGOLD™ is smaller than a single whole IgG molecule. In addition to a better penetration, the Fab' conjugates will have no interactions with Fc receptor-bearing membranes due to the absence of the Fc region in Fab' fragments.

Alexa Fluor® FLUORONANOGOLD™ is chromatographically purified through gel filtration columns. There are absolutely no aggregates or other molecular weight impurities. This is in sharp contrast to colloidal gold conjugates which usually are prepared by centrifugation to remove the largest aggregates, and frequently contain smaller aggregates.

Close to 1 NANOGOLD® label to 1 antibody or streptavidin makes this product distinct from the 0.2 - 10 variable stoichiometry of colloidal gold - antibody preparations.

Alexa Fluor® 647 FLUORONANOGOLD™ conjugates are used as a secondary antibody. They may also be used as a tertiary labeled antibody, or a custom Alexa Fluor®-FLUORONANOGOLD™ conjugate may be used as the primary antibody. If additional antibody incubation steps are used, rinse with buffer 3 (3 X 10 mins) after incubation.

Alexa Fluor® 647 FLUORONANOGOLD™ particles do not have affinity to proteins as do colloidal golds. This reduces background and false labeling.

Alexa Fluor® 647 FLUORONANOGOLD™ develops better with silver than do most colloidal gold giving it higher sensitivity. Both silver and gold enhancement can be used to make the immunolabeling useful for electron microscopy and light microscopy,

USING EM STAINS WITH Alexa Fluor® 647 FLUORONANOGOLD™

Because the 1.4 nm Alexa Fluor® 647 FLUORONANOGOLD™ particles are so small, over staining with OsO₄, uranyl acetate or lead citrate may tend to obscure direct visualization of individual NANOGOLD® particles. Four recommendations for improved visibility of Alexa Fluor® 647 FLUORONANOGOLD™ are:

1. Use of reduced amounts or concentrations of usual stains.
2. Use of lower atomic number stains such as NANOVAN™, a Vanadium based stain (Nanoprobes Catalog # 2011).⁵
3. Enhancement of Alexa Fluor® 647 FLUORONANOGOLD™ with silver developers, such as HQ SILVER™ (Nanoprobes Catalog#2012).
4. Enhancement of Alexa Fluor® 647 FLUORONANOGOLD™ with the gold developer, GOLDENHANCE™ (Nanoprobes Catalog#2113).

THIOL CAUTION

FLUORONANOGOLD[®] conjugates may be subject to loss of gold clusters (NANOGOLD[®]) upon exposure to thiols such as β -mercaptoethanol (BME) or dithiothreitol (DTT). Avoid use of thiol-based reducing agents. If a reducing environment is needed, reduce the protein, then purify from the thiol agent by column chromatography. Use non-metallic columns, and include 5 mM EDTA with the eluent, since trace metals catalyze thiol oxidation back to disulfides; most thiols do not reoxidize within several hours to several days following this procedure. Then use the Alexa Fluor[®] 647 FLUORONANOGOLD[™]. If a reducing agent is absolutely required, use a non-thiol agent, such as TCEP (tris(carboxyethyl) phosphine).

TEMPERATURE CAUTION

Although FLUORONANOGOLD[®] is stable under most conditions,⁶ labeled specimens or conjugates may not be stable above 80°C for long periods. Best results are obtained at room temperature or 4°C. It is best to use silver or gold enhancement before procedures requiring temperatures above 37°C, such as baking, or use low temperature embedding media (e.g., Lowicryl) if labeling before embedding.⁷

METHODS

Several publications describe the successful application of FLUORONANOGOLD[™] for light and electron microscopy. These provide additional protocols, details and applications that may be helpful in obtaining the best results (Refs. 1, 12-16).

FLUORESCENCE MICROSCOPY IMMUNOLABELING WITH Alexa Fluor[®] 647 FLUORONANOGOLD[™]

If aldehyde-containing reagents have been used for fixation, these should be quenched before labeling. This may be achieved by incubating the specimens for 5 minutes in 50 mM glycine solution in PBS (pH 7.4). Ammonium chloride (50 mM) or sodium borohydride (0.5 - 1 mg/ml) in PBS may be used instead of glycine.

The procedure below³ describes an example of the use of a Alexa Fluor[®] 647 FLUORONANOGOLD[™] conjugate as a secondary antibody probe. Dilutions of Alexa Fluor[®] 647 FluoroNanogold[™] will vary with different procedures, but a final concentration of 0.4 – 10 μ g/mL is advisable as a starting point for most applications; for simultaneous electron microscopy labeling, a compromise between the optimum concentrations for fluorescence and electron microscopy maybe necessary. Other protocols and techniques used with fluorescently-labeled antibodies may also be used with Alexa Fluor[®] 647 FLUORONANOGOLD[™] conjugates.

1. Fix cells in freshly-prepared 2 % formaldehyde in PBS for 15 mins at 20°C; alternatively, fix in 100 % methanol at -20°C for 3 minutes; if methanol fixation is used, skip to step 4.
2. Wash in PBS (3 x 10 mins).
3. Permeabilize in 0.2 % Triton X-100 plus 1 % normal serum (NS) from the host species of the Alexa Fluor[®] 647 FLUORONANOGOLD[™]-conjugated antibody in PBS at pH 7.4 for 5 minutes on ice.
4. Wash in PBS with 1 % NS (3 X 10 mins).
5. Incubate in the appropriate concentration of primary antibody for 1 hour at room temperature in a humidified chamber. If using 22 mm X 22 mm square cover slips, 30 μ L of diluted antibody is placed on the coverslip and the coverslip is inverted onto a glass slide. The slide is then placed in a humidified chamber which is incubated at room temperature.
6. Wash in PBS with 1 % NS + 5% non-fat dried milk (3 X 10 mins).
7. Incubate with Alexa Fluor[®] 647 FLUORONANOGOLD[™] reagent at a concentration of 0.4 – 10 μ g/mL (diluted in buffer containing 1% non-fat dried milk) for 1 hour in a humidified chamber at room temperature.
8. Wash in PBS (4 X 10 mins).
9. Mount coverslip with a drop of mounting medium. Observe as usual.

PBS Buffer:

20 mM phosphate
150 mM NaCl
pH 7.4

SUPERRESOLUTION MICROSCOPY WITH Alexa Fluor® 647 FLUORONANOGOLD™

A number of different superresolution methods are in use that have differing specimen preparation requirement. Since the requirements for each of these may be different, we do not provide a specific protocol. Instead, we recommend that you check with the manufacturer of your microscope system, or the literature in your field and area of application, for a current procedure and follow this.

General guidelines for stochastic optical reconstruction microscopy (STORM) are provided by Nikon, and may be found at:

<http://www.mvi-inc.com/wp-content/uploads/N-STORM+Protocol.pdf>

ELECTRON MICROSCOPY WITH Alexa Fluor® 647 FLUORONANOGOLD™

The procedures given in this section are complete immunolabeling procedures, and are also recommended for NANOGOLD® conjugates. If the specimen has already been labeled and observed by fluorescence microscopy, it requires only mounting, silver or gold enhancement (usually necessary) and staining according to your usual electron microscopy protocol before observation.

If aldehyde-containing reagents have been used for fixation, these should be quenched before labeling. This may be achieved by incubating the specimens for 5 minutes in 50 mM glycine solution in PBS (pH 7.4). Ammonium chloride (50 mM) or sodium borohydride (0.5 - 1 mg/mL) in PBS may be used instead of glycine.

Cells in Suspension

If the cells are already labeled, mount, stain and observe as usual. If a different specimen is to be used, the procedure below is recommended:

1. Optional fixing of cells: e.g., with glutaraldehyde (0.05 - 1% for 15 minutes) in PBS. Do not use Tris buffer since this contains an amine which reacts with glutaraldehyde.
2. Centrifuge cells (e.g. 1 mL at 10^7 cells/mL) at 300 X g, 5 minutes; discard supernatant; resuspend in 1 mL buffer. Repeat this washing (centrifugation and resuspension) 2 times.
3. Incubate cells with 0.02 M glycine in PBS (5 mins). Centrifuge, then resuspend cells in PBS-Milk buffer (specified below) or PBS containing 1 % BSA for 5 minutes.
4. Place 50 - 200 μ L of cells into Eppendorf tube and add 5 - 10 μ L of primary antibody (or antiserum). Incubate 30 minutes with occasional shaking (do not create bubbles which will denature proteins).
5. Wash cells using PBS-Milk as described in step 2 (2 X 5 mins). Resuspend in 1 ml PBS-Milk buffer.
6. Dilute Alexa Fluor® 647 FLUORONANOGOLD™ to a concentration of 0.4 – 10 μ g/mL in PBS-Milk buffer and add 30 μ L to cells; incubate for 30 minutes with occasional shaking.
7. Wash cells in PBS buffer as described in step 2 (2 X 5 mins).
8. Fix cells and antibodies using a final concentration of 1% glutaraldehyde in PBS for 15 minutes. Then remove fixative by washing with PBS buffer (3 X 5 mins).

PBS-Milk Buffer:

20 mM phosphate
150 mM NaCl
pH 7.4
1% Non-fat dried milk (final concentration)

Optional, may reduce background:

0.5 M NaCl
0.05% Tween 20
0.1% gelatin (high purity)

PBS Buffer:

20 mM phosphate
150 mM NaCl
pH 7.4

NEGATIVE STAINING

Negative staining may be used for electron microscopy of small structures or single molecules which are not embedded. Negative stain must be applied after the silver enhancement. NANOVAN™ negative stain is specially formulated for use with NANOGOLD® reagents;⁵ it is based on vanadium, which gives a lighter stain than uranium, lead or tungsten-based negative stains and allows easier visualization of Alexa Fluor® 647 FLUORONANOGOLD™ particles with little or no silver enhancement.

Thin Sections

Labeling with Alexa Fluor® 647 FLUORONANOGOLD™ may be performed before (the pre-embedding method)^{8,9} or after embedding and sectioning (the post-embedding method)^{8,9}. The procedures for both methods are described below.

Thin sections mounted on grids are floated on drops of solutions on parafilm or in well plates. Hydrophobic resins usually require pre-etching.

PROCEDURE FOR PRE-EMBEDDING METHOD:⁸

If specimen has already been labeled with Alexa Fluor® 647 FLUORONANOGOLD™, skip to step 9. If a fresh specimen is required for EM, the following procedure is recommended.

1. Float on a drop of water for 5 - 10 minutes.
2. Incubate cells with 1 % bovine serum albumin in PBS buffer at pH 7.4 for 5 minutes; this blocks non-specific protein binding sites and minimizes non-specific antibody binding.
3. Incubate with primary antibody, diluted at usual working concentration in PBS-Milk buffer or PBS containing 1 % BSA (1 hour or usual time. Buffer formulations are given below).
4. Rinse with PBS-Milk (3 X 1 min).
5. Incubate with Alexa Fluor® 647 FLUORONANOGOLD™ reagent diluted to a concentration of 0.4 – 10 µg/mL in PBS-Milk with 1 % normal serum from the same species as the Alexa Fluor® 647 FLUORONANOGOLD™ reagent, for 10 minutes to 1 hour at room temperature.
6. Rinse with PBS- Milk (3 X 1 min), then PBS (3 X 1 min).
7. Postfix with 1 % glutaraldehyde in PBS (10 mins).
8. Rinse in deionized water (2 X 5 min).
9. Perform silver or gold enhancement (e.g., HQ Silver™ or GoldEnhance™), as specified in those product directions.
10. Dehydrate and embed according to usual procedure.
11. Stain (uranyl acetate, lead citrate or other staining reagent) as usual before examination.

PROCEDURE FOR POST-EMBEDDING METHOD:⁸

1. Prepare sections on plastic or carbon-coated nickel grid. Float on a drop of water for 5 - 10 minutes.
2. Incubate with 1 % solution of bovine serum albumin in PBS buffer at pH 7.4 for 5 minutes to block non-specific protein binding sites.
3. Incubate with primary antibody, diluted at usual working concentration in PBS-Milk buffer or PBS containing 1 % BSA (1 hour or usual time. Buffer formulations are given below)
4. Rinse with PBS-Milk (3 X 1 min).
5. Incubate with Alexa Fluor® 647 FLUORONANOGOLD™ reagent diluted to a concentration of 0.4 – 10 µg/mL in PBS-Milk with 1 % normal serum from the same species as the Alexa Fluor® 647 FLUORONANOGOLD™ reagent, for 10 minutes to 1 hour at room temperature.
6. Rinse with PBS (3 X 1 min).
7. Postfix with 1 % glutaraldehyde in PBS at room temperature (3 mins).
8. Rinse in deionized water for (2 X 5 min).
9. If desired, contrast sections with uranyl acetate and/or lead citrate before examination.

Silver or gold enhancement may also be used to render the Alexa Fluor® 647 FLUORONANOGOLD™ particles more easily visible (see below); this is recommended if stains such as uranyl acetate or lead citrate are applied. Silver or gold enhancement should be completed before these stains are applied.

PBS-Milk Buffer:

PBS Buffer:

20 mM phosphate
150 mM NaCl
pH 7.4
1% Non-fat dried milk (final concentration)

20 mM phosphate
150 mM NaCl
pH 7.4

Optional, may reduce background:

0.5 M NaCl
0.05% Tween 20
0.1% gelatin (high purity)

SPECIAL CONSIDERATIONS FOR DIRECT VIEWING OF Alexa Fluor® 647 FLUORONANOGOLD™ IN THE ELECTRON MICROSCOPE

For most work, silver or gold enhancement is recommended to give a good signal in the electron microscope (see below). For particular applications, such as high resolution cryo-electron microscopy, visualization of the Alexa Fluor® 647 FLUORONANOGOLD™ directly may be desirable. Generally this requires very thin samples and precludes the use of other stains.

Alexa Fluor® 647 FLUORONANOGOLD™ provides a much improved resolution and smaller probe size over other colloidal gold antibody products. However, because NANOGOLD® is only 1.4 nm in diameter, it will not only be smaller, but will appear less intense than, for example, a 5 nm gold particle. With careful work, however, Alexa Fluor® 647 FLUORONANOGOLD™ may be seen directly through the binoculars of a standard EM even in 80 nm thin sections. However, achieving the high resolution necessary for this work may require new demands on your equipment and technique. Several suggestions follow:

1. Before you start a project with Alexa Fluor® 647 FLUORONANOGOLD™ it is helpful to see it so you know what to look for. Dilute the Alexa Fluor® 647 FLUORONANOGOLD™ stock 1:5 and apply 4 µL to a grid for 1 minute. Wick the drop and wash with deionized water 4 times.
2. View Alexa Fluor® 647 FLUORONANOGOLD™ at 100,000 X magnification with 10 X binoculars for a final magnification of 1,000,000 X. Turn the emission up full and adjust the condenser for maximum illumination.
3. The alignment of the microscope should be in order to give 0.3 nm resolution. Although the scope should be well aligned, you may be able to skip this step if you do step 4.
4. Objective stigmators must be optimally set at 100,000 X. Even if the rest of the microscope optics are not perfectly aligned, adjustment of the objective stigmators may compensate and give the required resolution. You may want to follow your local protocol for this alignment but since it is important, a brief protocol is given here:
 - a. At 100,000 X (1 X 10⁶ with binoculars), over focus, under focus, then set the objective lens to in focus. This is where there is the least amount of detail seen.
 - b. Adjust each objective stigmator to give the least amount of detail in the image.
 - c. Repeat steps a and b until the in focus image contains virtually no contrast, no wormy details, and gives a flat featureless image.
5. Now underfocus slightly, move to a fresh area, and you should see small black dots of 1.4 nm size. This is the NANOGOLD®. For the 1:5 dilution suggested, there should be about 5 to 10 gold spots on the small viewing screen used with the binoculars. Contrast and visibility of the gold clusters is best at 0.2 - 0.5 microns defocus, and is much worse at typical defocus values of 1.5 - 2.0 microns commonly used for protein molecular imaging.
6. In order to operate at high magnification with high beam current, thin carbon film over fenestrated holey film is recommended. Alternatively, thin carbon or 0.2% Formvar over a 1000 mesh grid is acceptable. Many plastic supports are unstable under these conditions of high magnification/high beam current and carbon is therefore preferred. Contrast is best using thinner films and thinner sections.

7. Once you have seen NANOGOLD[®] you may now be able to reduce the beam current and obtain better images on film. For direct viewing with the binoculars reduction in magnification from 100,000 X to 50,000 X makes the NANOGOLD[®] much more difficult to observe and not all of the gold particles are discernable. At 30,000 X (300,000 X with 10 X binoculars) NANOGOLD[®] particles are not visible. It is recommended to view at 100,000 X, with maximum beam current, align the objective stigmators, and then move to a fresh area, reduce the beam, and record on film.
8. If the demands of high resolution are too taxing or your sample has an interfering stain, or is thick, a very good result may be obtained using silver or gold enhancement to give particles easily seen at lower magnification.

SILVER ENHANCEMENT OF Alexa Fluor[®] 647 FLUORONANOGOLD[™] FOR EM

Alexa Fluor[®] 647 FLUORONANOGOLD[™] will nucleate silver deposition resulting in a dense particle 2-80 nm in size or larger depending on development time. It should be completed before any staining reagents such as osmium tetroxide, lead citrate or uranyl acetate are applied, since these will nucleate silver deposition in the same manner as gold and produce non-specific staining. Silver development is recommended for applications of Alexa Fluor[®] 647 FLUORONANOGOLD[™] in which stains are to be used, otherwise the Alexa Fluor[®] 647 FLUORONANOGOLD[™] particles may be difficult to visualize against the stain.

Our LI SILVER[™] silver enhancement system is convenient and not light sensitive, and suitable for all applications. Improved results in the EM may be obtained using HQ SILVER[™], which is formulated to give slower, more controllable particle growth and more uniform particle size distribution.¹⁰

Specimens must be thoroughly rinsed with deionized water before silver enhancement reagents are applied. This is because the buffers used for antibody incubations and washes contain chloride ions and other anions which form insoluble precipitates with silver. These are often light-sensitive and will give non-specific staining. To prepare the developer, warm the components to room temperature and mix equal amounts of the components immediately before use. Alexa Fluor[®] 647 FLUORONANOGOLD[™] will nucleate silver deposition resulting in a dense particle 2-20 nm in size or larger depending on development time. Use of nickel grids is sometimes preferred.

Fluorescence microscopy should be performed BEFORE silver enhancement. This is because the silver-enhanced gold particles can quench fluorescence.

The relevant procedure for immunolabeling should be followed. Silver enhancement is then performed as follows:

1. Rinse with deionized water (2 X 5 mins).
2. OPTIONAL (may reduce background)²: Wash several times with 0.02 M sodium citrate buffer, pH 7.0.
3. Float grid with specimen on freshly mixed developer for 1-12 minutes, or as directed in the instructions for the silver reagent. More or less time can be used to control particle size. A series of different development times should be tried, to find the optimum time for your experiment. With HQ Silver, a development time of 6 min. usually gives 15-40 nm round particles.
4. Rinse with deionized water (3 X 1 min).
5. Mount and stain as usual.

Fixing with osmium tetroxide may cause some loss of silver; if this is found to be a problem, slightly longer development times may be appropriate. Alternatively, use of 0.1 % osmium tetroxide instead of 1 % has been found to give similar levels of staining while greatly reducing etching of the silver particles.

NOTE: Treatment with osmium tetroxide followed by uranyl acetate staining can lead to loss of the silver enhanced NANOGOLD[®] particles. This may be prevented by gold toning:¹¹

1. After silver enhancement, wash thoroughly with deionized water.
2. 0.05 % gold chloride: 10 minutes at 4°C.
3. Wash with deionized water.
4. 0.5 % oxalic acid: 2 mins at room temperature.
5. 1 % sodium thiosulfate (freshly made) for 1 hour.
6. Wash thoroughly with deionized water and embed according to usual procedure.

GOLD ENHANCEMENT OF Alexa Fluor® 647 FLUORONANOGOLD™ FOR EM

The small 1.4 nm NANOGOLD® particles may alternatively be enhanced (grown to a larger size) for better visibility using GoldEnhance™, which catalytically deposits gold around the NANOGOLD®, making a larger solid gold particle. Gold enhancement may be preferable to silver enhancement in some cases due to the different properties of GoldEnhance™: a) Gold is chemically more stable and is not depleted by osmium or uranyl stains; b) Gold has higher backscattering and is useful for SEM; c) GoldEnhance™ is light insensitive – it can be used in normal room lighting, and development followed in the light microscope; d) GoldEnhance™ may be used with physiological buffers, such as ones containing chloride, which precipitates silver enhancers. GoldEnhance™ follows a similar procedure to silver enhancement, but for specific directions, see those that accompany GoldEnhance™.

REFERENCES

1. "Combined ALEXA-488 and Nanogold Antibody Probes;" Hainfeld, J. F.; Furuya, F. R.; Powell, R. D., and Liu, W.; *Microsc. Microanal.*, 8, (Suppl. 2: Proceedings) (*Proceedings of Microscopy and Microanalysis 2002*); Voekl, E.; Piston, D.; Gauvin, R.; Lockley, A. J.; Bailey, G. W., and McKernan, S., Eds.; Cambridge University Press, New York, NY, 2002, p. 1030CD (<http://www.nanoprob.com/applications/MSAFN02.html>).
2. Powell, R. D.; Hainfeld, J. F.; Halsey, C. M. R.; Joshi, V. N.; Hacker, G. W.; Hauser-Kronberger, C., and Takvorian, P. M.; *Microsc. Microanal.*, 5, (Suppl. 2: Proceedings); G. W. Bailey, W. G. Jerome, S. McKernan, J. F. Mansfield, and R. L. Price (Eds.); Springer-Verlag, New York, NY; 1999, 478-479 (<http://www.nanoprob.com/applications/MSAFN99.html>); Powell, R. D.; Halsey, C. M. R., and Hainfeld, J. F.; *Microsc. Res. Technique*, 42, 2-12 (1998); Powell, R.D., Halsey, C.M.R., Spector, D.L., Kaurin, S.L., McCann J., and Hainfeld, J.F.: *J. Histochem. Cytochem.* 45, 947-956 (1997).
3. Spector, D.L., and Smith, H. C.: Redistribution of U-snRNPs during mitosis. *Exp. Cell. Res.*, 163, 87-94 (1986).
4. Hainfeld, J. F., and Powell R. D.; *Cell Vision*, 4, 308-324 (1997); Furuya, F. R., and Hainfeld, J. F., *J. Histochem. Cytochem.*, 40, 177 (1992); Furuya, F. R., Hainfeld, J. F., and Powell, R. D., *Proc. 49th Ann. Mtg., Electron.Microsc. Soc. Amer.*; Bailey, G. W. (Ed.), San Francisco Press, San Francisco, CA; 1991, p. 284 (<http://www.nanoprob.com/applications/MSA92ng.html>).
5. Tracz, E., Dickson, D. W., Hainfeld, J. F., and Ksiezak-Reding, H. *Brain Res.*, 773, 33-44 (1997); Gregori, L., Hainfeld, J. F., Simon, M. N., and Goldgaber, D. Binding of amyloid beta protein to the 20S proteasome. *J. Biol. Chem.*, 272, 58-62 (1997); Hainfeld, J. F.; Safer, D.; Wall, J. S.; Simon, M. N.; Lin, B. J., and Powell, R. D.; *Proc. 52nd Ann. Mtg., Microsc. Soc. Amer.*; G. W. Bailey and Garratt-Reed, A. J., (Eds.); San Francisco Press, San Francisco, CA, 1994, p. 132 (<http://www.nanoprob.com/applications/MSANV.html>).
6. Hainfeld, J. F., and Furuya, F. R.; in *Immunogold-Silver Staining: Principles, Methods and Applications* (M. A. Hayat, Ed.), CRC Press, Boca raton, FL., 1995: pp. 71-96.
7. Krenács, T., and Krenács, L.; in *Immunogold-Silver Staining: Principles, Methods and Applications* (M. A. Hayat, Ed.), CRC Press, Boca raton, FL., 1995: pp. 57-69.
8. J. E. Beesley, in "Colloidal Gold: Principles, Methods and Applications," M. A. Hayat, ed., Academic Press, New York, 1989; Vol. 1, pp. 421-425.
9. Lujan, R.; Nusser, Z.; Roberts, J. D. B.; Shigemoto R.; Ohishi, H., and Somogyi, P.: *J. Chem. Neuroanat.*, 13, 219-241 (1997).
10. Humbel, B. M.; Sibon, O. C. M.; Stierhof, Y.-D., and Schwarz, H.: Ultra-small gold particles and silver enhancement as a detection system in immunolabeling and In Situ hybridization experiments; *J. Histochem. Cytochem.*, 43, 735-737 (1995).
11. Arai, R.; Geffard, M.; and Calas, A. Intensification of labeling of the immunogold silver staining method by gold toning. *Brain Res. Bull.*, 28, 343-5 (1992).
12. Robinson, J. M.; Takizawa, T.; Pombo, A., and Cook, P. R. Correlative fluorescence and electron microscopy on ultrathin cryosections: bridging the resolution gap. *J. Histochem. Cytochem.*, 49, 803-8 (2001).

13. Robinson, J. M.; Takizawa, T., and Vandre, D. D. Applications of gold cluster compounds in immunocytochemistry and correlative microscopy: comparison with colloidal gold. *J. Microsc.*, **199**, 163-79 (2000).
14. Takizawa, T., and Robinson, J. M. FluoroNanogold is a bifunctional immunoprobe for correlative fluorescence and electron microscopy. *J. Histochem. Cytochem.*, **48**, 481-6 (2000).
15. Takizawa, T., and Robinson, J. M. Analysis of antiphotobleaching reagents for use with FluoroNanogold in correlative microscopy. *J. Histochem. Cytochem.*, **48**, 433-6 (2000).
16. Takizawa, T., Suzuki, K., and Robinson, J. M. Correlative microscopy using FluoroNanogold on ultrathin cryosections. Proof of principle. *J. Histochem. Cytochem.*, **46**, 1097-102 (1998).

Technical Assistance Available.

For a complete list of references citing this product, please visit our world-wide-web site at

<http://www.nanoprob.com/references/Reffng.html>.